Table for section drying

OPERATING MANUAL

The undersigned authority legalizes the signature of the signing proxy(ies).

THE NETHERLANDS CHAMBER OF COMMERCE AND INDUSTRY

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Electronic Certifying Stamp

APOSTILLE

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1. Country: THE NETHERLANDS
This public document
2. has been signed by R.D.M. van den Dool
3. acting in the capacity of official of the
Chamber of Commerce and Industry and
4. bears the seal/stamp of aforesaid Chamber of
Commerce and Industry

Certified

5. in 's-GRAVENHAGE
6. on 1-12-2015
7. by the court registrar
8. no 2015-10526
9. Seal/stamp:
10. Signature:

E. van Houte

First we understand.

Then we innovate.

CONTENTS

1. INTRODUCTION1

 GENERAL INFORMATION1

 GENERAL DESCRIPTION1

 SAFETY INFORMATION1

 UNPACKING AND INSTALLATION.....1

 SCOPE OF DELIVERY.....1

2. OPERATION2

 2

 HEATING WORKING AREA2

 SETTING THE TEMPERATURE.....2

 CARE AND CLEANING.....3

3. SPECIFICATIONS4

1. Introduction

General information

Read this manual very carefully before using the equipment and always keep it handy. The equipment is not intended for use in explosive and/or flammable environments. Qualitative equipment operation can only be guaranteed when using original spare parts. The manufacturer is not liable for damage caused by incorrect operation.

General description

Heating stage is intended for drying paraffin sections placed on glass slides (this procedure may be required in histological, pathological, bacteriological, and chemical laboratories) after microtomy process (i.e., preparation of thin sections of the paraffin block using a specialized device - automated microtome).

Safety information

Heating stage complies with the European standards for electrical safety (EN 61010-1-1993). All metal parts of the stage which can come into contact with the user are grounded. Therefore, to ensure safe operation, turn on the stage into a grounded outlet. To enable electrical connection of the stage, use the supplied power cord. The equipment is not intended for use in potentially explosive atmospheres. The probability of an explosion can occur when using flammable substances (tools for anesthesia, skin care products and disinfectants). Maintenance should only be performed by trained service engineer. The manufacturer is not liable for damage caused by incorrect operation.

Unpacking and installation

Carefully unpack the stage and the accessories. When unpacking, do not use a knife or scissors to avoid damaging the equipment. Set the stage away from heat and direct sunlight, as this can lead to overheating. It is prohibited to work with the stage near the gas burners.

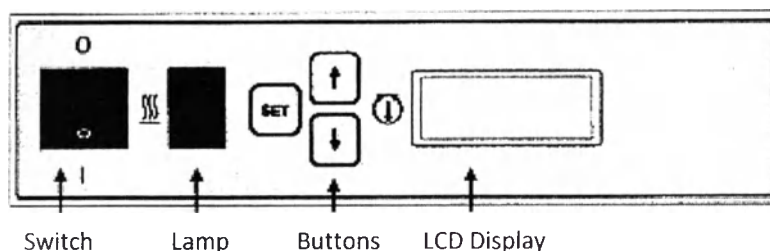
It is imperative that there is no dust and construction waste at the site of

installation/operation of the heating stage. Make sure that the mains voltage corresponds to the voltage indicated on the stage.

Scope of delivery

1. Table for section drying (article 1452)
2. Power cord
3. Operating manual.

2. Operation

Heating working area

WARNING: Before you start, check the set temperature: it should not exceed 60°C in order to avoid damage to the biopsy material.

1. Turn on the stage.
2. Press the switch. Lamp lights.
3. Set the desired temperature. The temperature is constantly displayed.
4. During the heating the lamp is green. It lights up when the required temperature is reached.

Setting the temperature

1. Press SET.
2. When the required temperature is saved, letter **P** appears on the screen.

3. By pressing “arrow up” and “arrow down”, you can increase or decrease the temperature by 1°C each time you press the button.

4. When you set the desired temperature, press SET to save it.
5. If during the temperature setting no button is pressed during 20 seconds, the stage turns off automatically.
6. The set temperature will be saved after the stage is turned off, as well as in case of power cuts.

Care and Cleaning

1. It is recommended to clean the stage every day.
2. Use only non-aggressive cleansers.
3. Cleaning powders can scratch the black stage coating.

NOTE: In case of damage of thermistor, **HLP** message appears on the screen and heating automatically stops. In this case, seek help from a service engineer.

3. Specifications

Dimensions	355 x 230 x 87 mm
Weight	3.9 kg
Voltage	230 V, 50 Hz
Power	460 V*A
Temperature range	30-70°C (adjusted by thermostat)
Temperature safety	> 90°C (protection against overheating)
Body material	PMS 427 C coated aluminum
Heating surface material	Black matte finish coated aluminum

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